



JAI Capture Interface 10 Installation Guide

Version: 10.0.0

Release Date: October 19, 2018

Installing the JAI Capture Interface


To acquire images with your imaging software, you must first install the appropriate capture interface. This is available from the Media Cybernetics support website at:

<http://www.mediacy.com/support/hardwaresupport> . These instructions cover the installation of the capture interface **only**.


You should also install the *eBUS SDK for JAI* camera software from *JAI*. This software may be downloaded from the JAI website .

To install the *JAI Capture Interface*, follow these steps.

1. Download the *JAI Capture Interface* from the *Media Cybernetics* support website and double-click on the installation application.
2. Click **Next** to begin installing the *JAI Capture Interface*.
You will see the *Welcome Screen*.

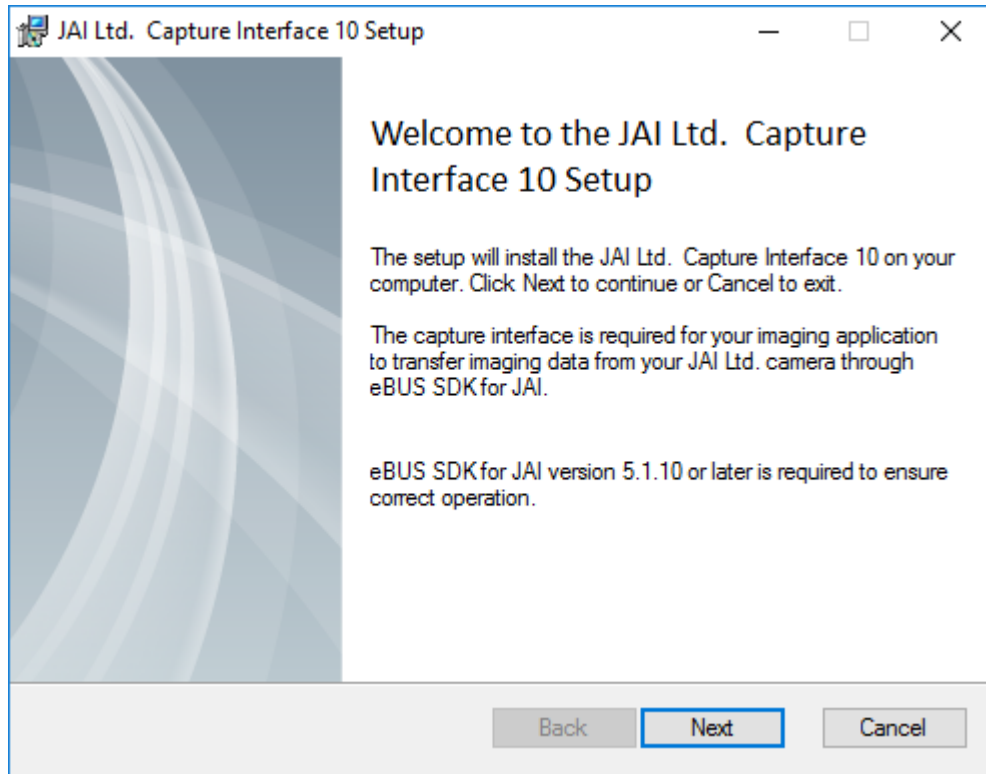


3. Click **Next** to begin installing the *JAI Capture Interface*.
You will see the *JAI Capture Interface* End-User License Agreement.

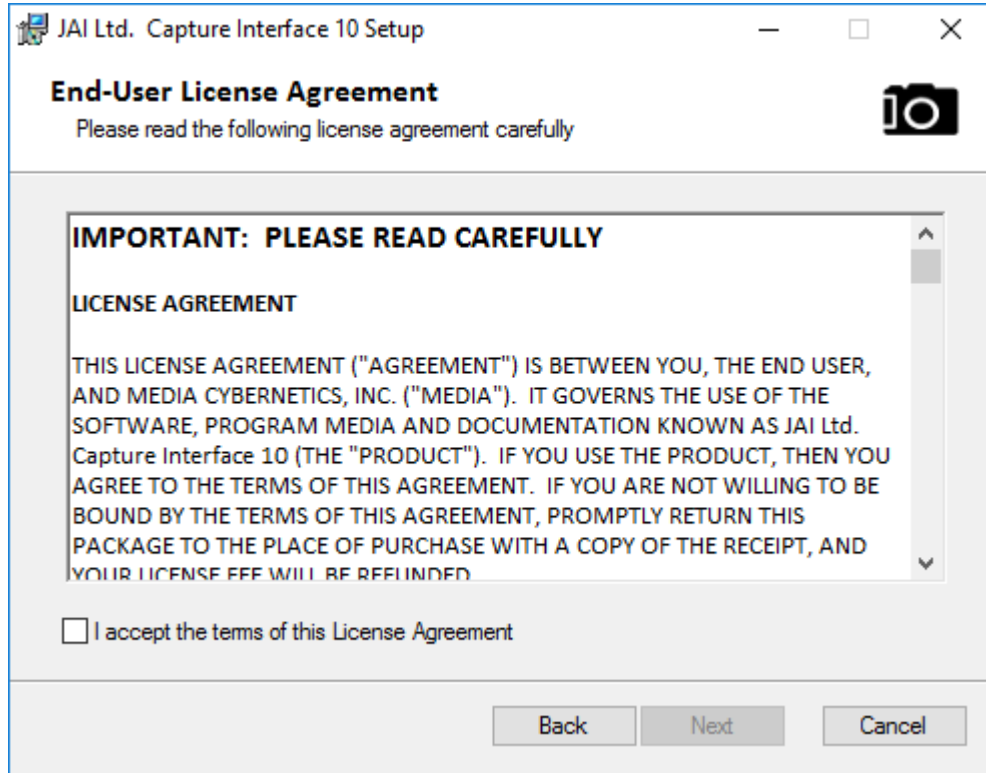


4. Check the box to accept the license agreement and click **Next**.


The *Select Target Application* screen appears.

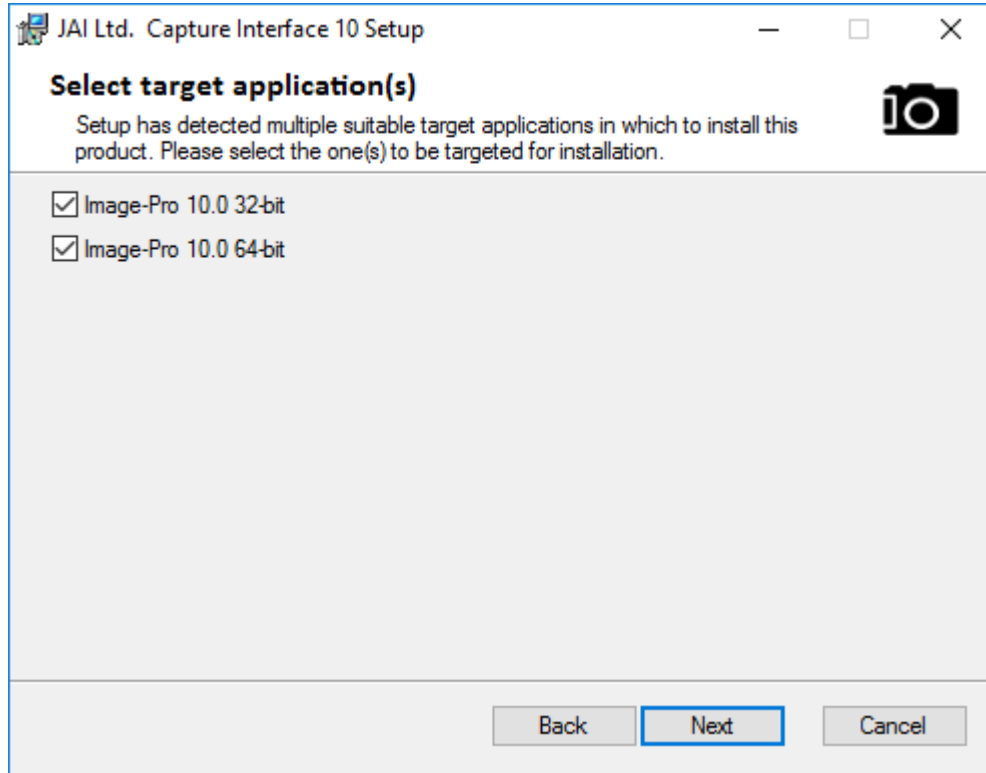

5. Check the box to indicate which software application you are using.
6. Click **Next** to continue.
The *Select Destination Folder* screen appears.

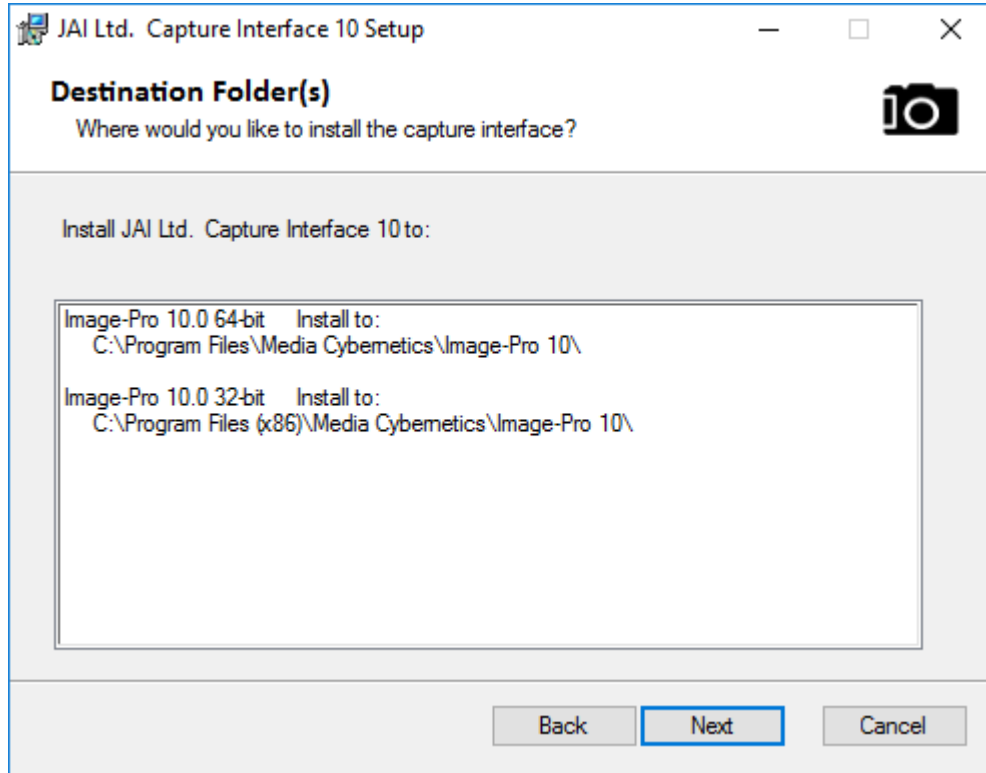


7. Indicate the program location, then click the **Next** button.
The *Ready to Install* dialog box appears.

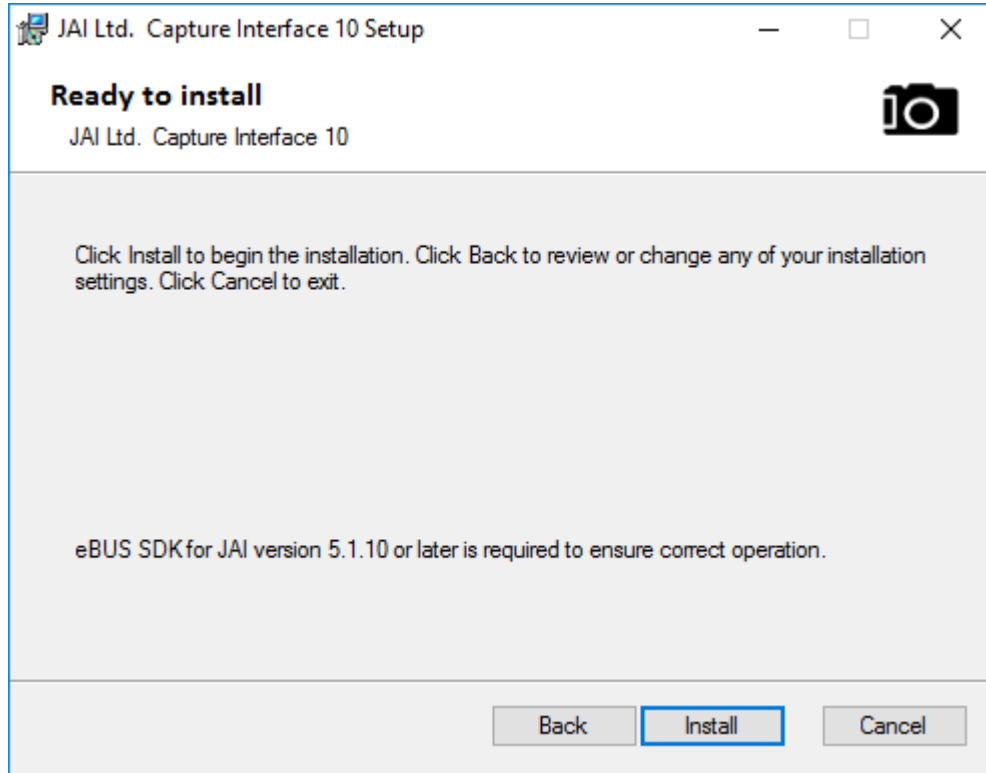

This screen indicates that your computer is ready to install the *JAI Capture Interface*.


8. Click **Install** to continue.

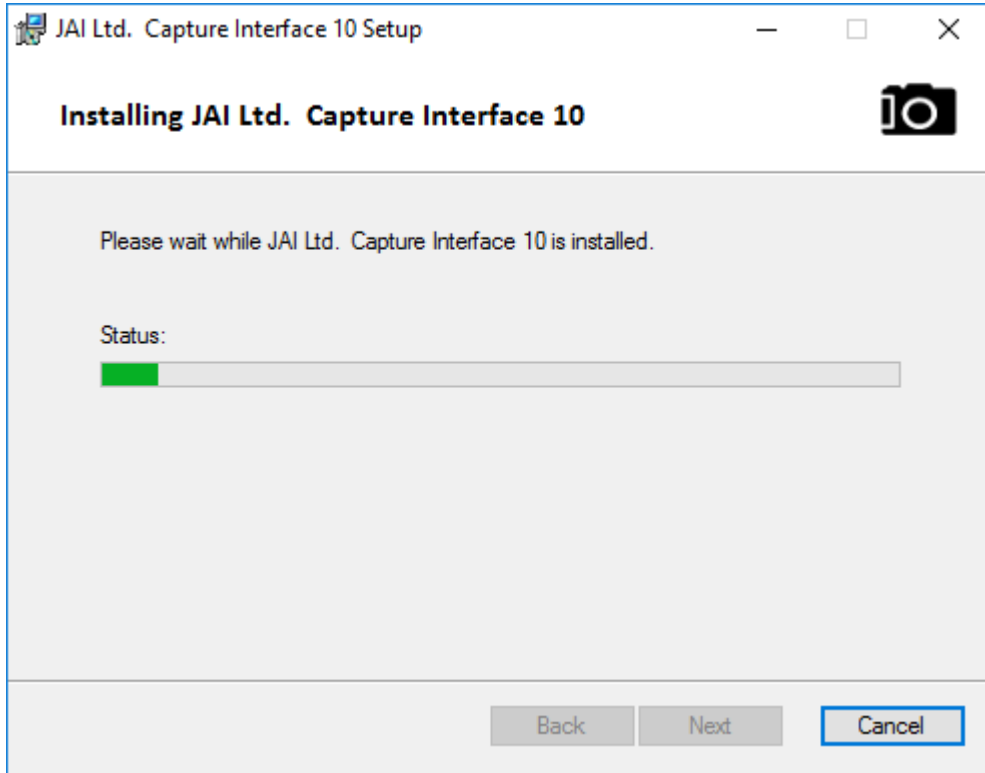

The *Installing* dialog box appears.

9. The setup program will begin copying and installing files. Click **Next** when the installation is complete.

The next page shows that the capture interface installation is complete.

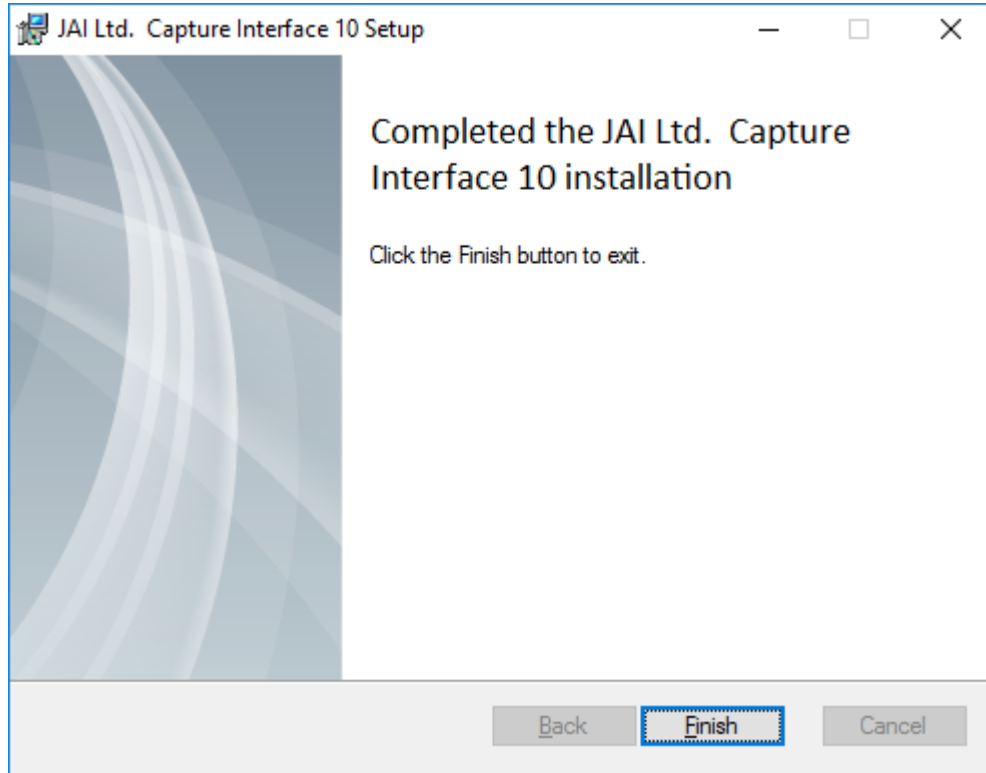

10. Click the **Finish** button.

You will see the following message:

