

## Installing the *QImaging PVCAM* Capture Interface

To acquire images with your imaging software, you must first install the appropriate capture interface from the Media Cybernetics support website:

<http://support.mediacy.com/drivers.asp>

These instructions cover the installation of the capture interface **only**.

You should also install the *PVCAM 2.9.11.3* driver provided by *QImaging*. The *QImaging* camera drivers can be downloaded from the *Media Cybernetics* website at


<http://www.mediacy.com/index.aspx?page=drivers> *QImg*

Recent improvements to the *QImaging* drivers may affect compatibility with older software products.


If your system is already configured to use your *QImaging* camera in another software application, please confirm compatibility with your software vendor prior to installing the latest version of the correct *QImaging* camera driver.

To install the *QImaging PVCAM* capture interface, follow these steps.

1. Click **Next** to begin installing the *QImaging* capture interface.

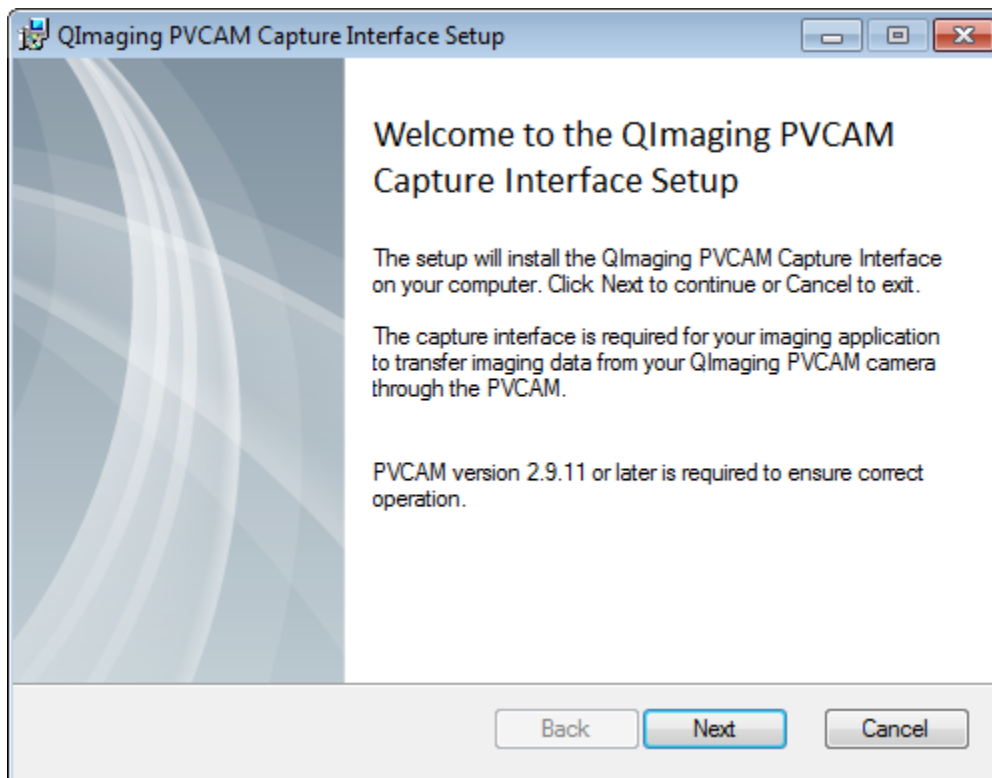


Next, you will see the *QImaging Capture Interface* End-User License Agreement.

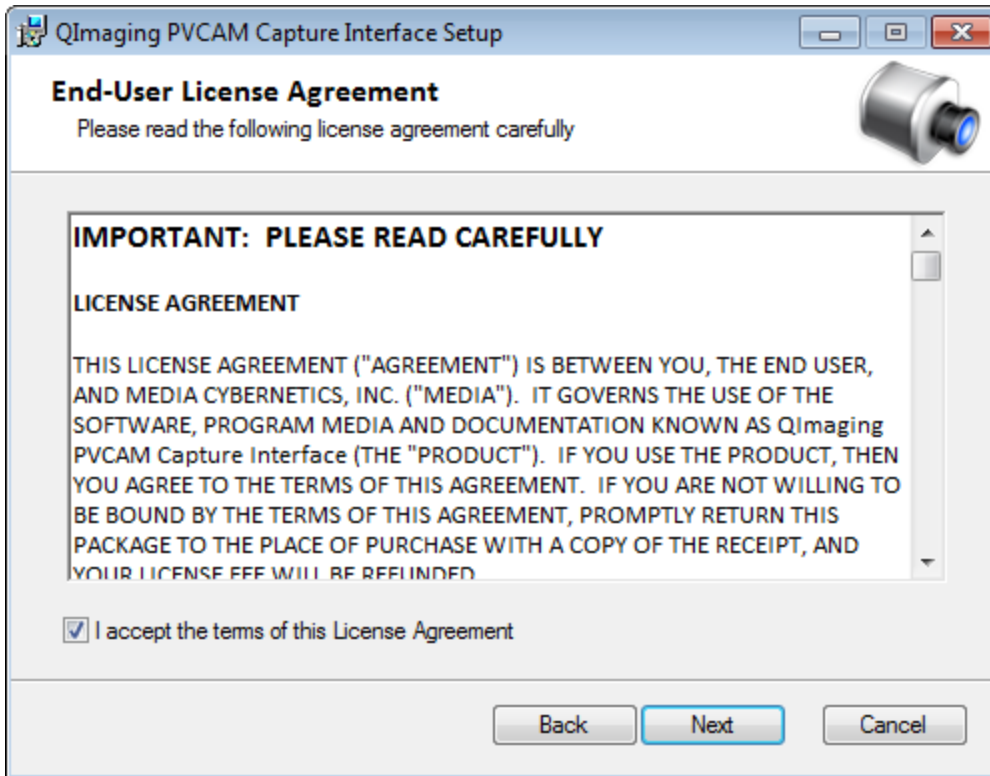


2. Check the box to accept the license agreement, and click **Next**.

The *Select Target Application* screen appears.

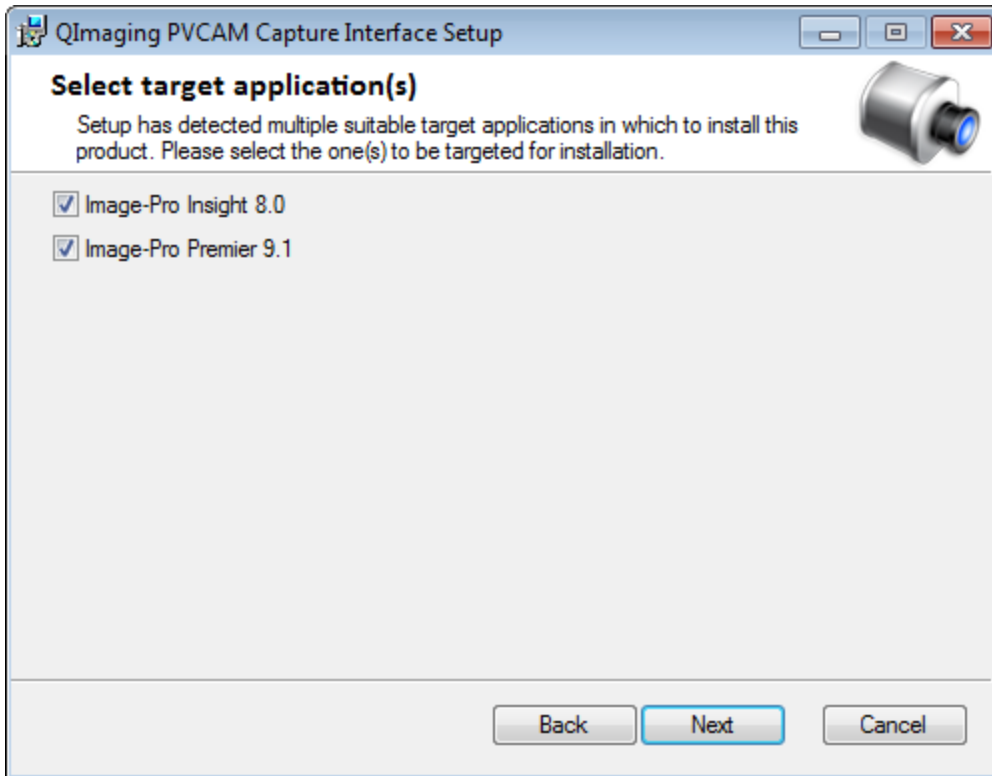


3. Check the box to indicate which software application you are using. Click **Next** to continue.

The *Select Destination Folder* screen appears.

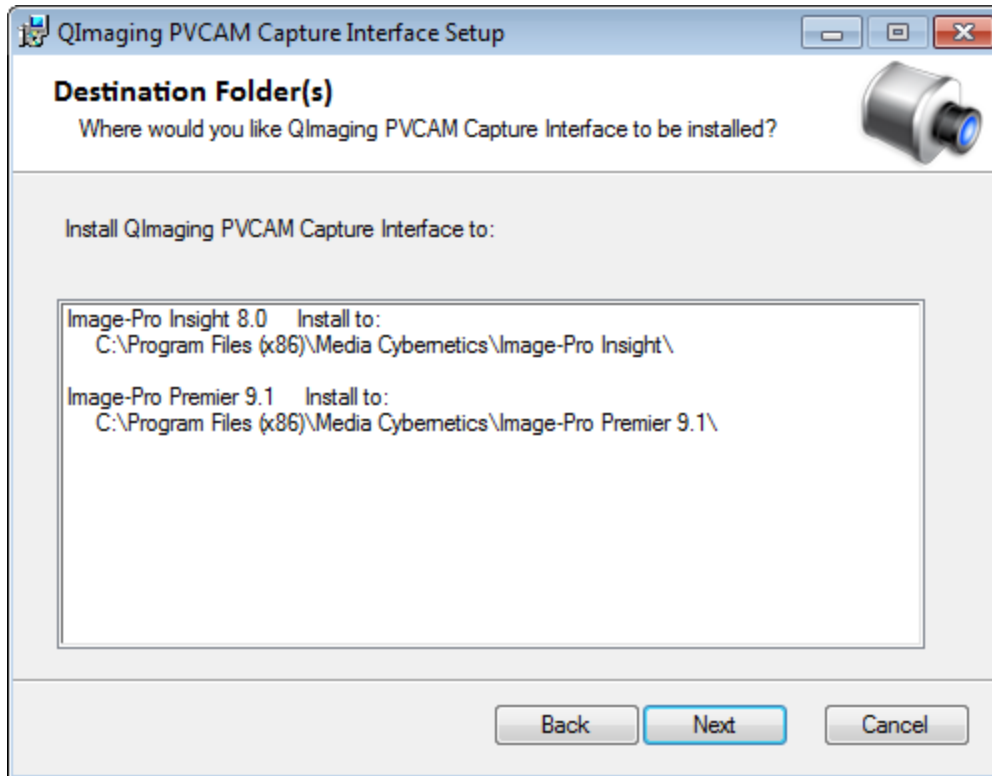

4. Indicate the program location, then click the **Next** button.


The *Ready to Install* dialog box appears.

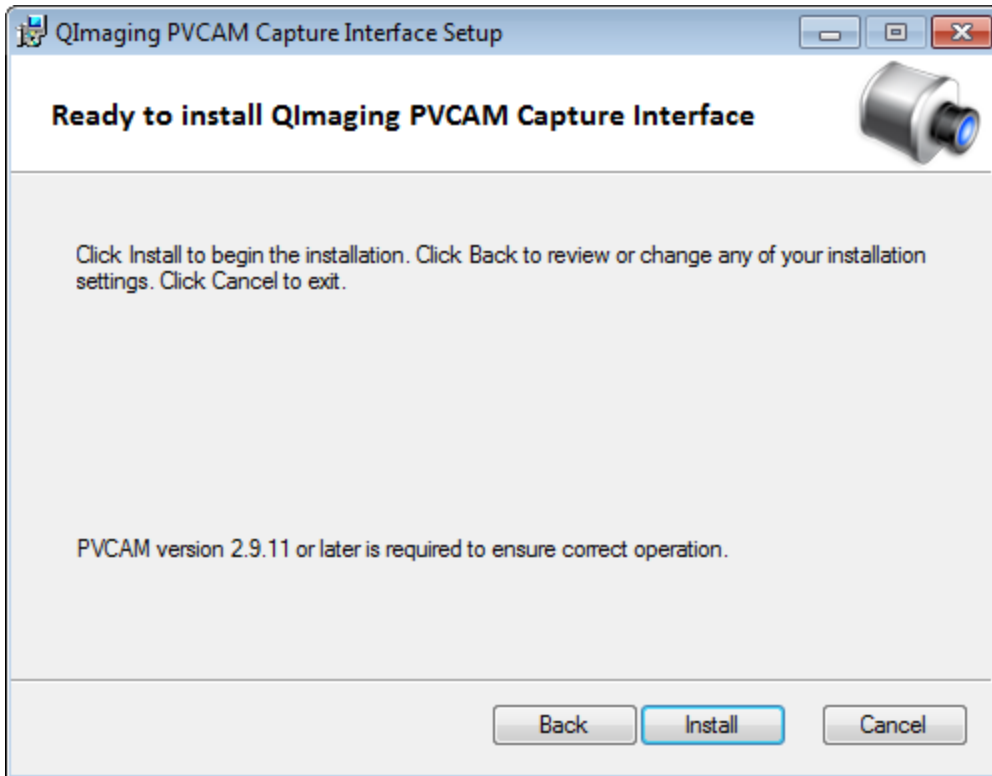

This screen indicates that your computer is ready to install the *QImaging PVCAM* capture interface.


5. Click **Next** to continue.

The *Installing* dialog box appears.

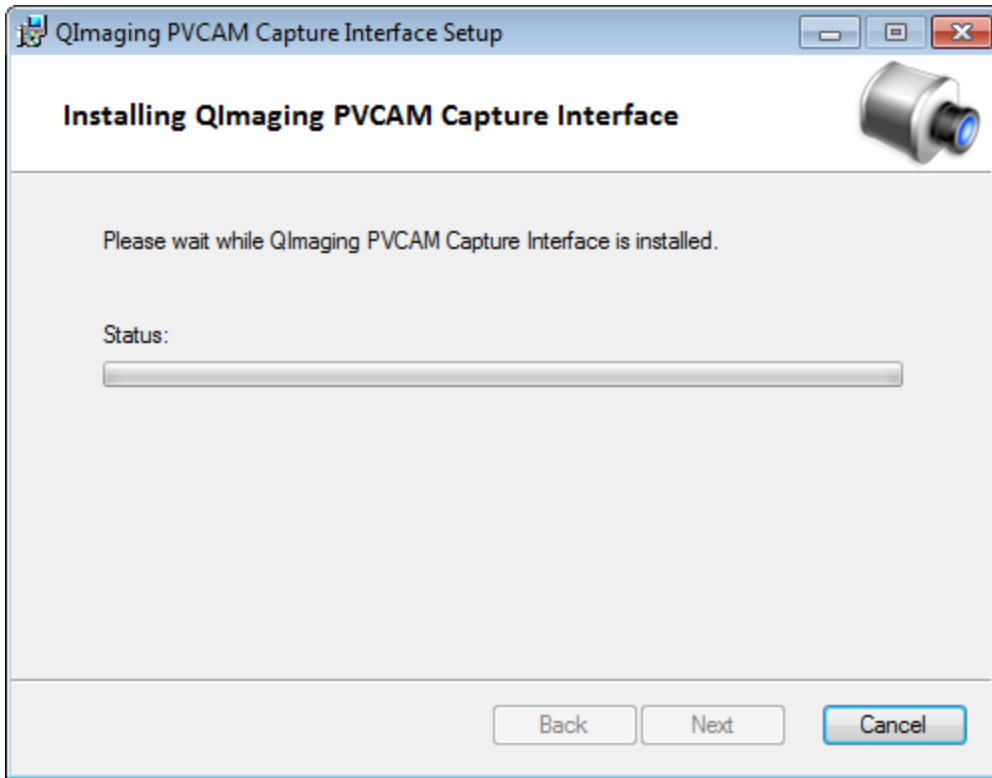


6. The setup program will begin copying and installing files. Click **Next** when the installation is complete.

The next page shows that the capture interface installation is complete.

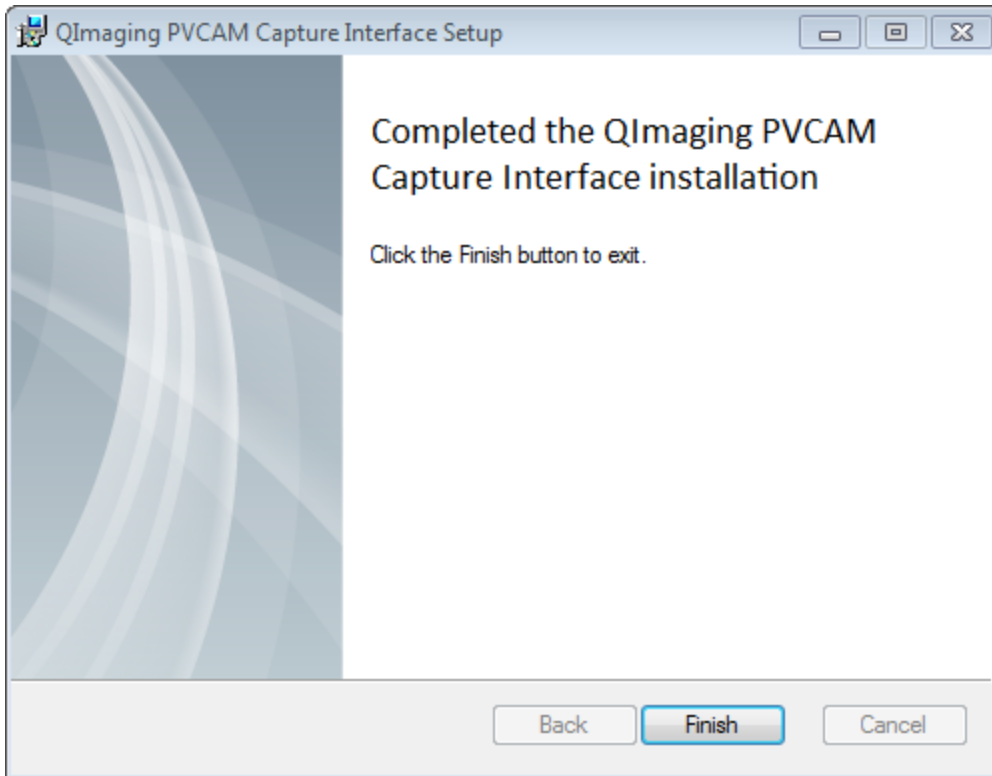

7. Click the **Finish** button.

You may see the following message:

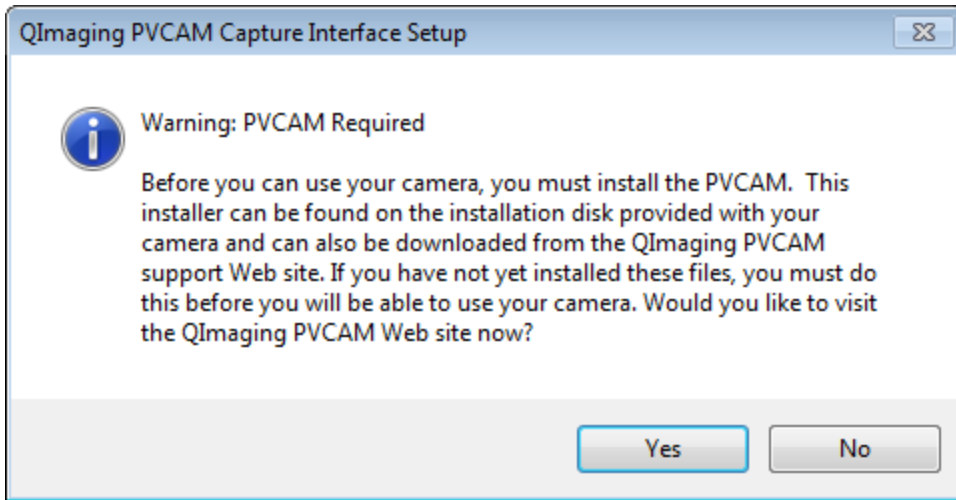

8. Click **YES** if you have not yet installed the PVCAM driver. You will be taken to the QImaging PVCAM website to complete the installation.